

# OPERATING MANUAL

## TABLE OF CONTENTS

**EchoTherm™  
PROGRAMMABLE  
DIGITAL HOT PLATE/STIRRER  
MODEL HS40/HS40A/HP40/HP40A  
Revised 9 January 2012**

**DOCUMENT NUMBER HS40-03**

TORREY PINES SCIENTIFIC, INC.  
2713 Loker Ave. West  
Carlsbad, CA 92010

TELEPHONE: (760)-930-9400

Toll Free: 866-573-9104

FAX: (760)-930-9480

E-MAIL: [info@torreypinesscientific.com](mailto:info@torreypinesscientific.com)

WEB SITE: [www.torreypinesscientific.com](http://www.torreypinesscientific.com)

	<u>PAGE</u>
I. INTRODUCTION	3
II. WARRANTY	3
III. RETURN OF ITEMS	3
IV. LABELS	3
V. CAUTIONS	4
VI. GENERAL DESCRIPTION	5
VII. FRONT AND REAR PANEL CONTROLS	6, 7
VIII. SET UP INSTRUCTIONS	8
IX. DISPLAY AND KEYBOARD DESCRIPTION	9, 10
X. SIMPLE OPERATIONS	11
XI. SIMPLE TEMPERATURE, RAMP RATE, STIRRER, TIMER, AND AUTO-OFF SETTINGS	11, 12
XII. PROGRAM OPERATIONS	13, 14
XIII. TEMPERATURE CALIBRATION	14, 15, 16
XIV. POWER INTERRUPTION PROTECTION	16
XV. RS232 INTERFACE	16, 17, 18
XVI. CLEANING, MAINTENANCE, AND CONSUMABLE PARTS	18, 19
XVII. ADDITIONAL SYMBOLS	19

## I. INTRODUCTION

Congratulations on your purchase of an EchoTherm™ Programmable Hot Plate or Stirring Hot Plate. Please read the instructions carefully to insure that you receive the maximum benefit from it. Also, be sure to fill out and return the enclosed warranty registration card.

## II. WARRANTY

Torrey Pines Scientific warrants this product to be free from defects in material and workmanship for a period of one year from the date of purchase. If repair or adjustment is necessary and has not been the result of abuse or misuse within the one year period, please return---freight prepaid--and correction of the defect will be made without charge.

Out-of-warranty products will be repaired on a charge basis.

## III. RETURN OF ITEMS

**Authorization must be obtained from our Customer Service Department before returning items for any reason.** When applying for authorization, please include data regarding the reason the items are to be returned. For your protection, items must be **carefully packed** to prevent damage in shipment and **insured** against possible damage or loss. Torrey Pines Scientific will not be responsible for damage resulting from careless or insufficient packing. A 15% restocking charge will be made on all unauthorized returns.

**NOTE:** Torrey Pines Scientific reserves the right to make improvements in design, construction, and appearance without notice.

## IV. LABELS

There are various labels on the body of this unit. Listed below are the labels and their meanings.

This symbol means: **Attention. The instruction manual is to be consulted for further information.** *Attention: Lire ce manuel pour des informations complémentaires.*



This symbol means: **Warning, Hot Surface.** *Attention, surface chaude.*



This symbol means: **Ground or earth connection.** *Connexion à la terre.*



This symbol means: **Dispose of in an environmentally correct manner.**

## V. CAUTIONS

### HEATER PLATE SURFACE

The Torrey Pines Scientific Programmable Hot Plates and Stirring Hot Plates are capable of temperatures in excess of 450°C at the plate surface. Touching the heated surface will cause severe burns. **USE EXTREME CAUTION AT ALL TIMES.** Never leave your hot plate accessible to others while it is hot. Although the unit is equipped with a **"HOT WARNING"** indicator do not rely on this alone. It is good practice to never touch the heater surface.

**NOTE: DO NOT INSTALL THIS UNIT CLOSER THAN SIX (6) INCHES (15.24 CM) TO A WALL OF COMBUSTIBLE MATERIAL. ALSO, THIS UNIT IS NOT FOR USE WITH FLAMMABLE SUBSTANCES.**

### TEMPERATURE PROBE

When attempting to control PROBE TEMPERATURE, plug in the temperature probe and place it in the sample AT ALL TIMES. If not placed into the sample and plugged into the rear of the hot plate, the unit will not be able to sense the temperature of the sample as heat is being applied. This results in driving the heater to its maximum and could result in ruining the sample.

Temperature probes vary in size and material. The most commonly used probe is the stainless steel 6-inch immersion probe (part number HS30-600). For other probes available in other lengths and materials, call the factory or consult your price list. All probes come with 3-foot (91.44 cm) interconnect cable and are 3/16th (47.6 mm) in diameter. Stainless steel probes work best with all but the most aggressive chemicals. Solid Teflon will work with the most aggressive chemicals, but only to about 260°C. Glass probes are available for use with aggressive chemicals at higher temperatures.

### ELECTRICAL

These hot plates and stirrers are made in models that operate at 100, 115, and 230 volts AC  $\pm 10\%$ . Be certain that your voltage matches the unit you receive. Check the nameplate on the bottom for the voltage setting on your unit. Be certain to use a three-wire, properly grounded AC input. Take the normal care and precaution one would use with any electrical appliance. Be careful to keep the AC line cord away from the hot plate. **Be certain to use a line cord of the same type and rating as the one supplied with this unit. Note: all fuses are fast blow. They are rated at 8 amps, 125 volts for ceramic topped 100 and 115 VAC units, and 4 amps, 250 volts for 230 VAC units. Cast aluminum topped units are 8 amps, 125 volts for 100 and 115 VAC units and 4 amps, 250 volts for 230 VAC units.**

*Bien vérifier que le cordon utilisé est du même type que celui livré avec l'unité. Note: tous les fusibles sont à action rapide. Lis sont à 8 A et 125 volts pour les unités à surface céramique de 100 et 115 volts et de 4A pour les unités de 250 et 230 volts. Pour les unités à boîtier en aluminium de 100 et 115 volts, l'ampérage est de 8A pour les fusibles. Pour les modèles en 250 et 230 volts, il est de 4A.*

## VI. GENERAL DESCRIPTION

The Torrey Pines Scientific Programmable Hot Plates and Stirring Hot Plates are general purpose digital hot plate/stirrer. They are available with a solid ceramic or cast aluminum heater tops. The top is 8" x 8" (20.32 cm x 20.32 cm) in size. All functions are settable from the digital front panel and display. These units when received display temperature in °C. They may be switched to read in °F. See page 6.

### HEATER

Either the plate surface temperature or the actual sample temperature may be set. A sensor in the plate is used to monitor surface temperature, or, alternately, a 100-ohm, 3-wire, platinum RTD temperature probe may be connected to the rear of the unit and inserted into the sample to measure sample temperature. When a temperature is set, power is applied to the heater to precisely control the temperature at the plate surface or at the sample, as directed by the user.

An optional "ramp value" may be entered into the unit which causes the temperature to approach the target value at a controlled rate of temperature change. This can be either an increase or decrease in temperature. Temperature is always displayed in degrees centigrade.

### TIMER

These units come with a count down timer that is settable and displays in hours, minutes and seconds. The timer can be set to 99:59:59 hours. When the timer is set it starts to count down to zero. At zero the timer will sound an audible alarm that will ring five times. In addition to the timer, these units have an AUTO-OFF function. When the AUTO-OFF function is activated the heater and stirrer will turn off at the end of any count down timer setting.

### ALARMS

In addition to the audible alarm associated with the count down timer, these units have an **over temperature alarm** that activates at 455°C (400°C on aluminum topped units). If, for any reason, the set temperature runs away and exceeds these limits, the unit will turn-on the over temperature alarm LED, sound an audible alarm, and shut the heater off. If or when the heater plate surface cools to below these limits, the heater will turn on again, and the alarm LED and audible alarm will turn off.


**NOTE: Failing to place the temperature probe in the solution and then setting a probe temperature will cause the heater to heat to over 455°C and set off the OVER TEMPERATURE ALARM. Also, the OVER TEMPERATURE ALARM may come on if a very large sample is placed on the unit and a probe temperature is set. This setting can cause the heater surface to go over 455°C to achieve the probe temperature set to heat the large sample.**

### STIRRER ( On stirring models only)

The stirrer is a motor driven magnet which revolves directly under the center of the heater plate. It is normal to stir a solution while heating. A stir bar is used in the solution and couples with the motor driven magnet and stirs the solution.

## VII. FRONT AND REAR PANEL CONTROLS

### FRONT PANEL

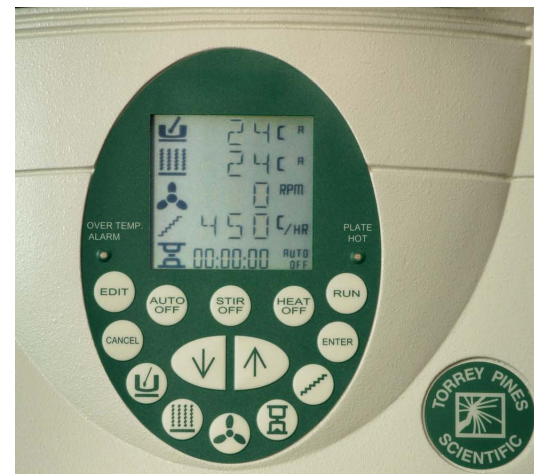


Shown above are the keyboard and display for the HS40 Programmable Digital Hot Plate/Stirrer. **(Note: The hot plate only models will not have the stirrer or stirrer off icons).**

The front panel has a tactile touch membrane keyboard with audible feedback. The keyboard is used to set all operating parameters. The display is a LCD type made up of five rows of numbers and several icons which function to display all parameters when the unit is running and during setting. When the unit is turned on, the display will come on and show the actual plate surface temperature. If a probe is used, it will display the probe temperature as well. The timer, stirrer and ramp functions will show zeros.

There are two LED's on the front panel as well. One is actuated when the plate surface goes over 50°C to remind the user that the plate surface is hot enough to cause burns. The second is for the over temperature alarm.

To change the unit from reading in °C to reading in °F or vice-versa hold the HEAT OFF button down for 3 seconds. The display will switch over. Degrees F is indicated when the "C" symbol next to the temperature reading for both plate and probe is missing. **The RAMP value is always set in °C/hour via the keypad or RS232 I/O port.**

## REAR PANEL

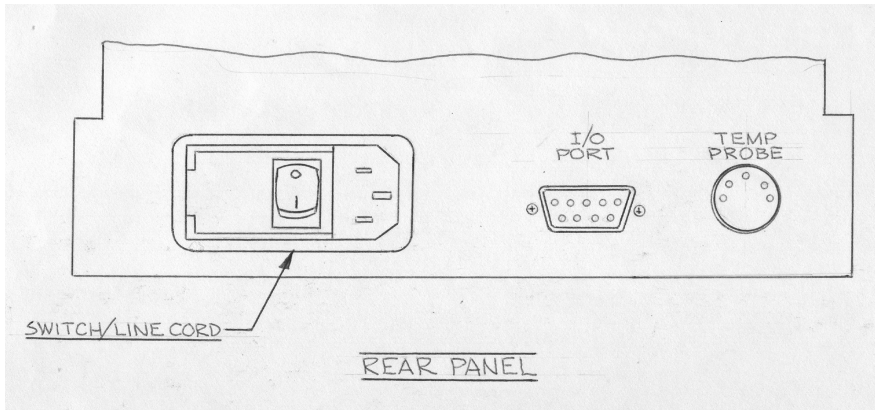

The AC power connector jack, fuse holder and switch are a module mounted on the right rear of the unit. The temperature probe jack is mounted to the right. Note the polarity on the probe jack. Be sure not to force this connection when plugging in the temperature probe.

The AC power jack is a three-prong, male plug combining the snap-in fuse holder and AC switch. The fuse types used appear on the label on the rear of the unit. Both neutral and high lines are fused.

**CAUTION: If the fuses blow repeatedly, contact your dealer or the manufacturer.**

**ATTENTION: Si les fusibles sautent de manière répétitive, contacter votre distributeur habituel ou le fabricant.**

The temperature sensor jack is a 3-pin DIN jack which is used with a 100-ohm at 0°C platinum RTD temperature probe.

## VIII. SET UP PARAMETERS AND INSTRUCTIONS

The HS40/HP40 Programmable units are a fairly simple instruments to install and use. Follow the instructions below.

### SET UP PARAMETERS

1. Ambient operating temperature range is from 5°C to 40°C.
2. Maximum altitude of operation should not exceed 2000 meters.
3. Maximum ambient operating relative humidity should not exceed 80% at 31°C decreasing linearly to 50% relative humidity at 40°C.

### SET UP INSTRUCTIONS

1. Place the unit on a level, dry bench or surface.
2. Plug the unit into a properly grounded, three-wire outlet of proper voltage.
3. Plug the temperature probe (if used) into the jack on the rear of the unit.
4. Place the sample on the heater plate and put the temperature probe (if used) into the sample container.
5. Turn on the unit by the switch on the rear panel. The unit will beep once and the display will light up. At this point the user can set or view any of the parameters of the unit.
6. Set either plate or probe target temperature, stirrer, timer, ramp rate, or program according to the instructions given in the following sections.

**Note: Do not use this equipment in any manner not specified by the manufacturer.**

**Note: Ne pas utiliser cet équipement si vous ne pouvez pas respecter les conditions d'utilisation spécifiées par le fabricant.**

### ENVIRONMENTAL INFORMATION

1. This unit is for installation category II.
2. This unit is rated pollution degree 2.

## IX. DISPLAY AND KEYBOARD DESCRIPTIONS

### DISPLAY

The display has five rows of numbers and icons. From the top they are as follows.

Probe Temperature: The icon to the left of the numbers represents a probe in a solution. This row of numbers shows the probe temperature. These numbers will always be followed with a "C" which denotes that the temperature reading is in degrees centigrade (unless in °F. In that case, the "C" will not appear next to the display). The small "T" and "A" following that tells the user that the numbers displayed are the actual temperature of the probe in the solution or the target temperature as set by the user. These two letters will toggle between TARGET(T) and ACTUAL(A), and the numbers in the display will change accordingly to show the progress of a set temperature toward its target. If no target temperature is set, then the display will always show the actual probe temperature.

Plate Temperature: The icon to the left of the second row of numbers represents the plate temperature at its surface as measured with a sensor in the heater plate. The row of numbers to the right of the icon shows the plate surface temperature. These numbers also are followed by a "C" which denotes that the temperature reading is in degrees centigrade (unless the display is reading in °F. In which case the "C" will not appear). The small "T" and "A" following that tells the user that the numbers displayed are the TARGET(T) or ACTUAL(A) plate surface temperatures. If no plate surface target temperature is set, the "T" and "A" will not toggle. However, the actual plate temperature is displayed even when a probe temperature has been set.

Stirrer (on stirring units only): The icon next in line down the display looks like a propeller and is the STIRRER icon. The row of numbers to the right of the icon displays the actual stirrer speed as set by the user. It is in RPM (revolutions per minute).

Ramp: The icon next in line down the display looks like a staircase and is the RAMP icon. The number value set in the display is always in degrees centigrade per hour. This tells the user that the target temperature as set for the probe or plate is moving from its starting point to the target temperature at a fixed rate in degrees centigrade per hour.

Timer: The next icon down the display is the hourglass icon that represents the timer function. The numbers that can be set and displayed next to the icon are in hours, minutes, and seconds. The timer can be set to 99:59:59 maximum. The timer will count down to zero from its setting and sound an audible alarm and then start to count up letting the user know how long it has been since the timer counted down to zero. The words "AUTO-OFF" next to the numbers are visible only when the AUTO-OFF function is actuated. When this function is actuated, the timer will count down to zero, sound the audible alarm, and turn the heater and stirrer off. In addition, the Timer numbers are used in writing a program, and in displaying program steps and cycles as the program progresses. More on writing and running programs later in the manual.

### KEYBOARD

Up/Down Arrows: The UP and DOWN arrows in the center of the keyboard are used to set values for the probe temperature, plate temperature, stirrer, ramp, timer, and progressing within the steps of a program when writing or editing the program. Pushing and holding the UP arrow will cause the display parameter as selected to scroll up in value. Pushing and holding the DOWN arrow will cause the display parameter as selected to go down in value. In addition, the DOWN arrow is used as "cancel" to turn off the beep when the timer counts down.

Probe Temperature Icon: The probe temperature icon is the one that matches the icon to the left of the top line of the display. Touching and releasing this icon selects the UP and DOWN arrows to set the PROBE TEMPERATURE.

Plate Temperature Icon: The plate temperature icon is the one that matches the second line down from the top in the display. Touching and releasing this icon selects the UP and DOWN arrows to set the PLATE TEMPERATURE.

Stirrer Icon (on stirring units only): The stirrer icon matches the third line down of the display. Touching and releasing this icon selects the UP and DOWN arrows to set the STIRRER value.

Ramp Icon: The ramp icon matches the fourth line down from the top of the display. Touching and releasing this icon selects the UP and DOWN arrows to set the ramp value.

Timer Icon: The timer icon is the one that matches the last line at the bottom of the display. Touching and releasing this icon selects the UP and DOWN arrows to set the TIMER value.

Auto-Off Key: Touching and releasing this key activates the AUTO-OFF function. When activated, this will shut off the heater and stirrer when the timer counts down to zero. This function can be turned on or off by touching the AUTO-OFF button at any time during a timed event. When activated, the words "AUTO-OFF" will be illuminated on the display.

Heat Off Key: Touching and releasing the HEAT OFF key will turn off the heater. The HEAT OFF key also is used to change the unit temperature display from °C to °F and vice-versa. Holding the key down for 3 seconds will cause the display to change.

Stir Off Key (on stirring units only): Touching the STIR OFF key turns off the stirrer.

Edit Key: The EDIT key is used to enter the program memory to either write or edit a program.

Cancel Key: The Cancel key is used to cancel a program step while writing or editing a program.

Run Key: The Run key is used to start running a program in memory.

Enter Key: The Enter Key is used to enter a value into a program.

## X. SIMPLE OPERATION

These are fully programmable units. With them user has the ability to write and store in memory ten programs of as many as ten steps each for instant recall and later use. Also, the unit can be used to set a single temperature, temperature ramp, stirring speed (on stirring units) and timer function without going into the memory of the unit. The memory should be used to store complex routines where all the parameters of the sample can be changed automatically to other parameters without user attention.

### XI. SIMPLE TEMPERATURE, RAMP RATE, STIRRER, TIMER, AND AUTO-OFF SETTINGS

#### TEMPERATURE

The heater may be set to control either the plate surface or the sample itself. This is done by entering a target temperature from the front panel keyboard. The control electronics in the hot plate will automatically apply power to the heater plate to reach the desired temperature.

The user may enter either a target probe temperature or a target plate temperature. Only one target temperature is allowed at one time. Setting a target probe temperature will erase any target plate temperature previously set and vice versa. Target temperatures may be set anywhere in the range from ambient to 450°C (842°F) or 400°C (752°F). The unit comes with the display reading in degrees C. To change to degrees F depress and hold the HEAT OFF button until the display changes. To change back to degrees C repeat the process. When in degrees F the C symbol next to the temperature reading will not display.

Setting Solution Temperature: To set a solution temperature, touch and release the solution temperature icon. Next, touch the UP arrow until the display (top row) shows the temperature wanted. If the target is overshoot while setting, push the DOWN arrow until the correct value is in the display. The unit will now turn the heater on and drive the solution to the target temperature. Note that the solution display will toggle between the TARGET (T) and the ACTUAL (A) temperatures. When the target temperature is reached, the "T" and "A" will alternate but the numerical value will remain the same.

**CAUTION: Be certain the temperature probe is in the solution and plugged into the rear panel jack when a solution temperature target has been entered. Failure to do so could damage your sample because the hot plate will drive to maximum while seeking a temperature it cannot find.**

**PRÉCAUTION D'EMPLOI: Bien vérifier que la sonde de température plonge dans le liquide et qu'elle est connectée dans la fiche située à l'arrière de l'appareil lorsque vous sélectionnez une température. L'échantillon risque d'être endommagé si la sonde n'est pas connectée car la plaque va monter au maximum de sa température en recherchant une température qui n'aura pas été.**

Setting Plate Temperature: To set a plate temperature, touch and release the plate temperature icon. Next, press the UP arrow until the display (second row) shows the temperature wanted. If the target is overshoot while setting, push the DOWN arrow until the correct value is in the display. The unit will now turn on the heater and drive the plate to the target temperature selected. Note that the display will toggle between the TARGET (T) and ACTUAL (A) temperature. Again, when the target temperature is reached, the "T" and "A" will alternate but the numerical value displayed will remain the same. **Note: When a heater value is turned off, the value is saved so that when the heater is next used that previously set value is displayed and set. To change that value use the Up or Down arrows.**

Setting Ramp: **NOTE: IF A RAMP VALUE IS TO BE USED, THE RAMP VALUE MUST ALWAYS BE SET BEFORE SETTING A TARGET TEMPERATURE.** To set a ramp value, touch and release the ramp icon. Press the UP arrow until the display (fourth row) shows the value desired. The heater will now go from its current value to the target temperature, either plate or solution, at the rate just instructed. The ramp value is always in degrees centigrade per hour even if the temperature display has been set to read in degrees F. The RAMP rate is settable from 1°C/Hour to 450°C/Hour. Note that the unit will heat fastest when **not** using a ramp rate setting, even 450°C/Hour. The ramp rate should only be used to slow the heating rate.

It is valuable to note that the ramp can be set to go up to a target temperature or down from a higher temperature to a lower one. However, the unit can never ramp down any faster than the sample can cool naturally. When ramping down, the heater is used to slow a cooling rate only.

Setting Stirrer (on stirring units): To set a stirrer speed, touch and release the stirrer icon. Press the UP arrow until the display (third row) shows the value desired. The numbers shown are in revolutions per minute (rpm). The stirrer will now go to the value instructed.

Setting Timer: To set a timer value, touch and release the timer icon (hourglass). That icon will now flash in the display. Note that the seconds digits in the display will flash. Seconds can now be set if wanted by pressing the UP arrow. To set minutes, press the timer icon again. Note that the minute digits will now flash. Minutes can now be set by touching the UP arrow. To set hours, press the timer icon again and the hours digits will flash. Hours can now be set by touching the UP arrow. If only seconds or minutes are set, continue to press the timer icon until it scrolls past hours. Touching the timer icon again will start the timer. When the Timer counts down to zero it will alarm 5 times and start to count up. The up counter lets the user know how long it has been since the Timer counted to zero. To stop the timer from counting up, touch the DOWN arrow.

Setting Auto-Off: To set the auto-off function, touch and release the AUTO-OFF button. The display will illuminate the words "AUTO-OFF" to the right of the timer telling the user that the auto-off function is set. To turn off the auto-off function, touch the AUTO-OFF button again. The words, "AUTO-OFF" will no longer be illuminated in the display. Remember, when AUTO-OFF is on, the unit will automatically turn off the heater and stirrer when the timer counts down to zero.

## XII. PROGRAMMABLE OPERATION

The HS/HP40 Series have the ability to store ten programs in memory for instant recall and use at any time. Each program has ten steps maximum where one step is a temperature, stirring speed, ramp function (if wanted), and time. At the end of each program there is a cycle step that can be selected which will repeat the program up to 98 times. Setting the cycles to 99 makes the program repeat infinitely. These programs are stored in CMOS and cannot be lost by a power failure or by turning the unit off. However, if power fails during a program run, when power returns, the unit will have lost its instructions and will be sitting with the same information on the display as when the unit was first turned on. To stop and program while it is running touch CANCEL.

### Some rules for programming:

**To erase all programs in memory, depress and hold the EDIT key for about 5 seconds. The display will be all blank with the exception of 4 small lines at the bottom of the display. When the EDIT key is released, after a second or so, the display will return to the normal screen and the programs will be erased. This erases ALL programs in memory. To change or erase one program the user must go into the program and remove each step.**

**The timer can be used in two ways in a program. The timer can be used with a ramp function or without a ramp function. When the timer is used with a ramp function, the timer does not start to count down until the target temperature is reached. When the timer is used without a ramp function, the timer starts to count down as soon as the program starts. If the user wants a timed event to start after the target temperature is reached, but does not want to wait for a ramped approach to the target temperature to take place, set the ramp to 450. In almost all cases, this ramp rate is faster than the unit can approach the target temperature, and it will appear as though no ramp rate is being used, and the timer will start to count down after the target temperature is reached.**

### RUNNING A PROGRAM STORED IN MEMORY

1. Press RUN. The display will show PR 01 at the bottom of the display. This is program 1.
2. If program 1 is the choice to be run, next press ENTER. Program 1 will now start to run. If program 1 is not the choice use the UP ARROW to scroll to the program of choice. When the program is chosen, press ENTER and the program stored in that site will start to run. As the program starts to run, as the unit progresses through the steps in the program, the audible alarm will sound once for step 1, twice for step 2, three times for step 3, etc.

## HOW TO WRITE A PROGRAM

**Remember, a program step is a temperature, stirring speed (if wanted), temperature ramp (if wanted), and a timer setting. Do not advance to the next step until the current one is complete.**

- 1 Press EDIT
- 2 Select the program site (1 through 10) using the UP and DOWN ARROWS.
- 3 Touch ENTER. The display will show 01 01, for program one step one.
- 4 Enter the ramp value if wanted (remember how the timer works with and without a ramp function) and touch ENTER. Then enter a temperature and ENTER, and stirring speed and ENTER, and the time function for that step and ENTER. This step is now finished.
- 5 Press the UP ARROW to go to step 2. The display will show 01 02 for program one step two. Now enter the values for ramp, temperature, stirrer and timer being certain to hit ENTER after each. Step 2 is now complete.
- 6 Repeat the steps above until the program is written.

**Note: If a mistake is made while entering a program step, just hit CANCEL on that step and make the new entry. To go back to a prior program step, hit the DOWN ARROW. Likewise, to go forward to another program step, hit the UP ARROW.**

- 7 After the last program step data has been entered, press ENTER twice. The display will now show CYC 01. If only one cycle of the program is needed, hit ENTER. If more than one cycle of the program is wanted, use the UP ARROW to go to the number of cycles wanted. For example, twenty cycles would appear on the display as CYC 20. Then press ENTER again and the program is written and stored for use.
- 8 To run a program, press RUN, the display will show PR 01. To run program 1, hit ENTER. To select and run a different program, use the UP ARROW to scroll to the program to be run, and press ENTER.

## XIII. TEMPERATURE CALIBRATION

The HS/HP40 Series are designed for accuracy. The temperature calibration designed into the unit is made to hold for very long periods of time. When calibrated in the factory, it is expected that the unit will meet the most demanding customers' requirements. However, our standards for temperature measurement may not be the same as the users. Therefore, these units have been designed to be calibrated in the field by the user. Follow the instructions below if calibration is wanted or needed.

**Note:** The units come calibrated from the factory, however, changes to any calibration point or points can still be made in the field. The procedure is listed below. To calibrate the unit the user is advised to remove all previously entered calibration data. That is done by depressing and holding the AUTO-OFF key until the second beep. Then depress and hold either the PLATE or PROBE key (depending on which is to be erased and calibrated) until the second beep. This will erase all calibration data previously stored in memory in either the PLATE or PROBE calibration. Follow the instructions below if calibration is wanted or needed.

#### PROBE CALIBRATION

Probe calibration is performed by using an accessory calibration kit which precisely simulates fixed temperature points. The kit has two "dummy" probes which, when plugged into the probe jack, represent the value of resistance equal to the temperature shown on the dummy probe. The dummy probes represent temperatures of 25°C and 400°C. These are the probe calibration points. The kit is available from the manufacturer (HS30-700).

To calibrate the probe temperature follow these steps.

1. Insert the 25°C dummy probe in the probe jack at the rear of the unit. Depress and hold the AUTO-OFF key for 3 seconds followed by the PROBE key and the display will flash. Use the UP or DOWN arrow as needed to set the display to read 25°C. The press and release the PROBE key again. When the display stops flashing it will read 25°C and the low calibration point has been set.
2. Remove the 25°C dummy probe and replace it with the 400°C dummy probe. Depress and hold the AUTO-OFF key for 3 seconds followed by the PROBE key. The display will flash. Use the UP or DOWN arrow to make the probe display read 400°C. When the display stops flashing it will read 400°C and the high calibration point has been set.
3. Remove the dummy probe. The probe calibration is now complete.

#### PLATE CALIBRATION

The plate calibration does not affect the probe calibration and vice versa. If you are using the probe temperature control loop only, it is not necessary to calibrate the plate loop. The plate loop has been factory calibrated and is not likely to change.

To calibrate the plate, one needs an accurate temperature meter, preferably digital, and a surface temperature probe. This is a very difficult measurement to make, and not all surface temperature probes will do the job well. If you need help, please contact the factory. The Plate and Probe calibration kit #HS30-800 is available from the factory and will contain the temperature meter, probe, and dummy probes for the probe calibration.

The calibration procedure requires that calibration starts at the low temperatures and goes high. It is time consuming. Be sure before making an adjustment that the plate temperature has stabilized at the point being calibrated. Follow the procedure below.

1. Start at ambient. Using your temperature meter, read the plate surface temperature in the center of the plate. Push and hold the AUTO-OFF key for 3 seconds followed by the PLATE key. The plate display will flash. Use the UP or DOWN arrows to make the display agree with the measured plate temperature on your meter. Then press and release the PLATE key again. When the display stops flashing the low calibration point is set.
2. Set the PLATE temperature to 400°C. The plate will start to heat. Give the unit ample time to reach and stabilize at 400°C. When the temperature is stable, measure the plate temperature with your meter. Depress and hold the AUTO-OFF key for three seconds until the plate display flashes. Use the UP or DOWN arrow as needed to make the plate temperature display read what is measured on your meter. When the display stops flashing the high temperature calibration point is set.

The plate surface calibration is now complete.

#### **XIV. POWER INTERRUPTION PROTECTION**

The unit is provided with power interruption protection. If AC power is interrupted while running a temperature, the unit will go back to the set temperature when power returns. To let the user know that a power interruption has occurred, the temperature display will be flashing between TARGET(T) and ACTUAL(A) faster than usual by only dwelling for one second on each. Under normal operating conditions, the display will alternate between (T) and (A) dwelling on (A) for three seconds and (T) for one second. To restore the display to normal, touch the DOWN arrow.

#### **XV. RS232 INTERFACE**

The RS232 is available through the 9-pin D-subminiature connector on the rear of the unit. It operates at 2400 Baud, 8 Data Bits, 1 Stop Bit, No Parity. No handshake hardware or software is necessary. It will work well on a Windows terminal program per the settings listed here. All communications settings and queries are done using ASCII characters with carriage return as the terminating character (CR<sup>1</sup>). **The more complete serial command set is at the rear of this manual.**

Any serial port on your PC can be used. The pin outs follow standards for RS232C. On the connector, pin 5 is ground, pin 3 is RX and pin 2 is TX. We do not provide software for this application, however, several users have found that using Windows Terminal Program, and Visual Basic works well, but any programming language can be used.



**Note: To comply with CE and to avoid possible EMI radiation from the RS232 cable, you should use a shielded cable.**

<u>Query</u>	<u>Command</u>
Request current plate temperature	a
Request current probe temperature	b
Request current timer	c
Request ramping value	d
Request plate target temperature	e
Request target probe temperature	f
Request stirrer target speed (rpm)	g

For example: To request current plate temperature one needs to send two characters

a + CR

and the unit responds xxx + CR where xxx is the present temperature of the heater plate.

NOTE: If there is no target plate temperature set into the unit it will send "---" + CR.

<u>Set</u>	<u>Command</u>
Plate target temperature	A
Probe target temperature	B
Timer	C
Ramping	D
Stirrer speed (rpm)	E
Stirrer Off	F
Heater Off	G
Toggle Auto On/Off	H

For example: To set the following command string do the following:

To set the plate temperature target to 123°C

A 123 + CR

To set the timer

C 12:34:5600 + CR

To set stirrer speed to 500 rpm

E 500 + CR

Note: If the command string is valid, the unit will respond back with the message "Command OK" + CR. If the command is incorrect the message will be "Command Failed" + CR.

## **XVI. CLEANING, MAINTENANCE, AND CONSUMABLE PARTS**

### **Cleaning**

This unit is subject to splashes and spills during normal use. Be sure to clean all spills quickly. Wipe spills with a soft cloth or paper towel. If a cleaning solution is necessary, use a mild soap or detergent solution and a soft cloth. Be sure not to use solvents. They could damage the paint, labels or display window on the unit. A mild abrasive can be used to clean the ceramic glass heater surface. Do not use abrasives on the cast aluminum heater top. It will scratch the surface.

**CAUTION: Do not attempt to clean the heater surface when it is hot. Burns could occur.**

***PRÉCAUTION D'EMPLOI: Ne pas nettoyer la surface de la plaque lorsqu'elle est chaude pour éviter toute brûlure accidentelle.***

### **Maintenance**

There is no ongoing maintenance program needed with this unit other than the normal care and cleaning as instructed above, and a simple inspection done whenever the unit is to be used. This simple inspection should include:

- 1- Checking that the AC line cord is not frayed or burned.
- 2- Checking that the unit is not dirty to a point where proper performance is impaired. This is especially important relative to the membrane switch and LCD window.
- 3- Being certain to store the unit properly, when not in use, in an area that will not have items placed on top of the unit, and covered in a way that will keep dirt and other foreign bodies out of the unit.

### **Spare and Consumable Parts**

There are very few spare or consumable parts necessary. A simple list with part number is below.

<u>Part Number</u>	<u>Description</u>
HS30-710	Replacement Keyboard Covers
730-0001	Power Cord, US
730-0006	Power Cord, German (European)
730-0008	Power Cord, UK
730-0004	Power Cord, Italian
730-0005	Power Cord, Australian
380-0003	Fuse, 8 amp, 125 volt
380-0004	Fuse, 10 amp, 125 volt
380-0005	Fuse, 4 amp, 250 volt

### **XVII. ADDITIONAL SYMBOLS**

The following are additional symbols found on the labels of the instrument.

<u>Symbol</u>	<u>Description</u>
V	Voltage
~	Alternating Current
A	Current
Hz	Frequency
W	Power
F	Fast Acting Fuse
T	Time Delay Fuse
I	Mains On
O	Mains Off

# Serial Interface Specification

---

for the

## Torrey Pines Scientific Models HP30, HS30 and HS40 Series Hotplates

Revision A

Torrey Pines Scientific  
2713 Loker Ave. West  
Carlsbad, CA 92010

Tel 760-930-9400  
Fax 760-930-9480

[www.torreyпинессcientific.com](http://www.torreyпинессcientific.com)

### About this Specification

This document describes the operation of the serial interface port on the Torrey Pines Scientific Model xx Hotplate. The serial interface is used to read and set values, such as the temperature or stirrer speed, from a computer.

The specification is intended for expert software programmers who wish to write low-level drivers for the hotplate.

### Hardware Description

The serial interface port uses standard RS-232 voltage levels and utilizes three wires: one for data from the computer to the hotplate, one for data from the hotplate to the computer, and ground.

Signal	9-Pin Serial Port
Data from Hotplate to Computer	Pin 2
Data from Computer to Hotplate	Pin 3
Ground	Pin 5

*Pinout and Cable Wiring*

A standard serial cable may be used to connect the hotplate to the computer.

The transmission of the data bytes on the serial interface is at 2400 baud, with no parity, 8 data bits, and one stop bit.

### Character Strings

Data is sent to and from the hotplate in ASCII character strings. Data exchange is always initiated by the computer. The computer sends a character string to the hotplate and the hotplate sends a string back to the computer in response. A response is always generated by the hotplate when it receives a carriage return.

Command String:     <Command> <Data> <cr>

Response String:     <Data> <cr>

The command defines whether and what type of data is required in both the command and response strings.

Each string is terminated by a carriage return character. Characters may be sent to the hotplate with any amount of delay between characters. The characters are accumulated into a buffer and held until the carriage return character is received. Once the hotplate receives the carriage

return character in a command string, it responds immediately. Character strings are sent from the hotplate with no delay between the characters.

*Note - no handshaking signals are used. Synchronization of the computer and the hotplate is accomplished through the command/response sequence only.*

*Note - be sure not to send the hotplate extra characters, such as linefeeds, or subsequent commands may be misinterpreted.*

## Commands

The specific commands and the associated data flow is shown in the table below:

Command Code	Command	Data Sent	Data Returned	Data Format
a	Get Plate Temperature	none	Plate Temperature	TEMP
b	Get Probe Temperature	none	Probe Temperature	TEMP
c	Get Timer	none	Timer	TIMER
d	Get Ramp	none	Ramp Value	TEMP
e	Get Plate Target	none	Plate Target	TEMP
f	Get Probe Target	none	Probe Target	TEMP
g	Get Stirrer Speed	none	Stirrer Speed	RPM
A	Set Plate Target	Plate Target	Result	TEMP
B	Set Probe Target	Probe Target	Result	TEMP
C	Set Timer	Timer	Result	TIMER
D	Set Ramp	Ramp Value	Result	TEMP
E	Set Stirrer Speed	Stirrer Speed	Result	RPM
F	Set Stirrer Off	none	Result	--
G	Set Heater Off	none	Result	--
H	Toggle Auto Off Mode	none	Result	--

Command Set

## Data Formats

The various data types are defined below.

### TEMP

Temperature in °C. (For the ramp value, this is °C per hour.)

### RPM

Stirrer speed in RPM.

### TIMER

Timer in hours, minutes, and seconds (hh:mm:ss).

## Result String

The result string returned from the hotplate for the "set" commands is one of three strings: "Command OK" if the operation is successful, "Command Failed" if the command cannot be executed, or "Invalid Command" if the command code is not in the list above.

The command will fail if the data is out of the acceptable range or in an invalid format.

## Examples

- To read the current plate temperature, send:

a<cr>

A typical response may be:

123<cr>

meaning that the plate is at 123°C.

- To set the stirrer speed to 300 r.p.m., send:

E300<cr>

The response will be:

Command OK<cr>

and the stirrer will turn on and go to 300 r.p.m.

- To set the timer to 5 minutes and 12 seconds, send:

C00:05:12<cr>

The response will be:

Command OK<cr>

and the timer will start running at 5:12.

